Evap-Trap 120 ml EX
Instructions

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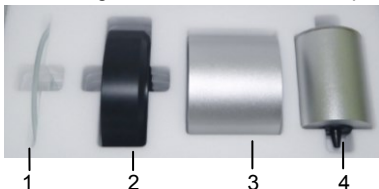
KIT CONTENT

The following items are included in the package:

1. Glass Plate Cover
2. Evaporation Trap (45 ml)
3. Wind Ring
4. Reservoir (120 ml)

I. ASSEMBLY INSTRUCTIONS

Follow the steps below. Refer to the figures when performing these steps.

Note: It is highly recommended to remove all draft-shield doors to allow for easy access while pipetting.

- 1** Remove the glass shelf.

- 2** Remove the side doors.

- 3** Remove the weighing pan.

- 4** Remove the balance wind ring.

- 5** Install the reservoir into the load cell's mounting cone.

- 6** Install the wind ring, making sure the side with the indentation is facing up.

- 7** Install the evaporation trap and then add water to the evaporation trap's trench, making sure that the water level is below the lower trench wall so it does not spill over.

Note: Max. 45ml, recommended water amount is 20~35ml

Add water here

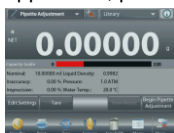
- 8** Install the glass plate cover. After all parts have been installed properly, it should look like this:

II. PIPETTE ADJUSTMENT APPLICATION

After the assembly is finished, tare the balance by pressing the **Tare** button. Enter the Pipette Adjustment Application on the Explorer balance and follow the screen instructions. For detailed instructions on how to use this application, please refer to the Explorer instruction manual.

III. USING THE PIPETTE

Dispense the water through the hole in the top glass cover plate. When dispensing the water with the pipette, make sure to hold the pipette at an angle, as shown below, for best results. For more information on how to use the pipette, please refer to the pipette instruction manual:

<http://dmx.ohaus.com/WorkArea/showcontent.aspx?id=4294968610>

- [Ohaus AO pipette instruction manual \(EN\)](#)

IV. TECHNICAL DATA

Model	Pipette Adjustment Kit
Oder Number	30252145
Applications	Pipette Adjustment
Trap Capacity	Maximum 45 ml
Reservoir/Container	120 ml
Minimum Dispensed Volume	> 100 μ l
Shipping Weight	1.1 kg with carrying case
Shipping Dimensions (W x L x H)	1.3 in x 0.6 in x 1.3 in / 33 mm x 16 mm x 33 mm

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P/N 30278712C © 2016 OHAUS Corporation, all rights reserved/ todos los derechos reservados / tous droits réservés / Alle Rechte vorbehalten / tutti i diritti riservati.
OHAUS Corporation